



DESIGN. INTEGRITY. RELIABILITY.

RODENT SLING FRAME & COVER: INSTRUCTIONS FOR USE



Figure 1: Rodent Sling Frame & Cover



Figure 2: Rodent Sling Frame & Suit

FRAME:

The rodent sling frame is routinely supplied with a one-size adjustable canvas and mesh cover. This is shown in the picture above (Figure 1). The design allows for animals to be quickly located in the sling and for easy adjustment and fit between individual animals.

SLING SUIT:

The rodent sling suit (purchased separately) is best described as a full body rat jacket with canvas extensions through which the supports of the frame are passed (Figure 2).

The body portion of the suit is made from our lycra/spandex rat jacket fabric and closes by means of a Velcro insert. This provides for a quick, secure fit and adjustment between individual animals.

The extensions are made from canvas fabric with a sleeve on each side through which the bars of the frame are passed.

The overall dimensions make the suit compatible with the sling frame and the jacket portion of the suit corresponds to a full body rat jacket.

It is recommended that the animals be acclimated to the suit in advance of any procedures.

Should you have any questions or need more information please contact our Customer Service Department toll free in North America at 1-877-425-3604 or send an e-mail to info@lomir.com.

LOMIR Biomedical Inc.

CANADA:
USA

95 Huot, Notre-Dame-de-l'Île-Perrot, QC Canada J7V 7M4 / Tel: (514) 425-3604 / Fax: (514) 425-3605 / Toll Free: 1-877-425-3604

444 East Main Street, Malone, NY, USA 12953 / Tel: (518) 483-7697 / Fax: (518) 483-8195 / Toll Free: 1-877-425-3604

Email: info@lomir.com / Website: www.lomir.com /