


The logo for Lomir Biomedical Inc. features the word "Lomir" in a white, sans-serif font. The letter "o" is stylized with a white arc that curves over the top of the letter and extends to the left, ending in a small dot. The background is a dark blue gradient with a pattern of lighter blue circles of varying sizes.

BIOMEDICAL INC.

CANADA | USA | EUROPE | LOMIR.COM

Restraint Slings Features and Uses

RESTRAINT SLINGS

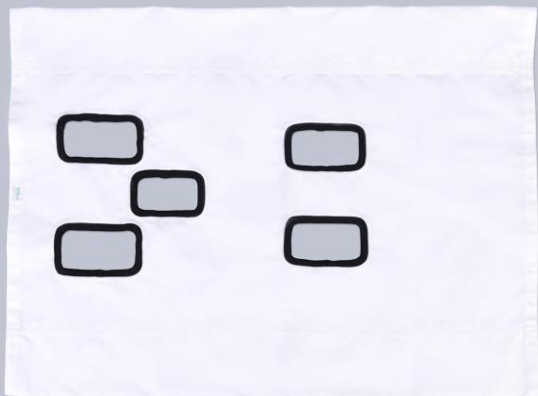

- Comfortable and secure restraint is essential when investigations call for animals to be held in devices for any length of time
- LOMIR manufactures sling restraints for various applications and species

LARGE ANIMAL FRAMES

FRAME



- The frame is constructed from 304 stainless steel 1" x 1" x .65" square tube. The complete frame is welded which provides a rigid construction. Dimensions for a standard frame are 24" wide x 36" long x 39" high. The frames are mounted on 5" diameter castors.
- The three foot length frame is suitable for dogs, mini-pigs, juvenile farm pigs and small ruminants.
- The four foot length frame being for large dogs, swine and farm animals.
- The deep "V" at the front of the frame allows easy access for instrumentation, sampling, administration and examinations.



LARGE ANIMAL FRAMES

SLING COVER



- The versatility of the frame is increased simply by having other sling covers on hand; covers are available for specific weight ranges and for a variety of species.
- The sling cover is made from a high quality coated canvas for durability. All openings and are edged in cotton jersey for comfort. Additional fabrics may be used for slings and custom features, such as openings for data recording can be added.

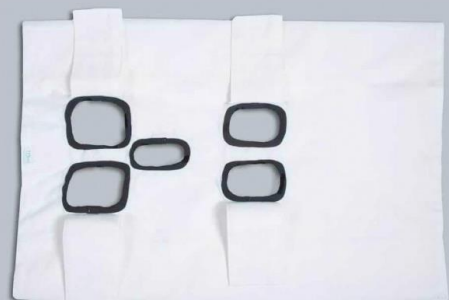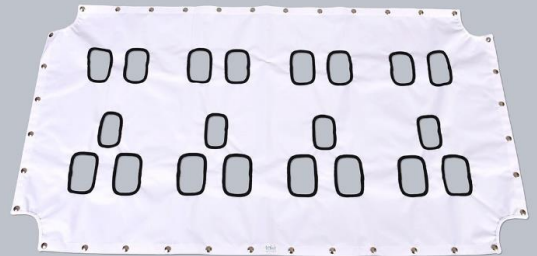

LARGE ANIMAL FRAMES

FEATURES


- The frame features two locking castors. Foot operation holds frame in place.

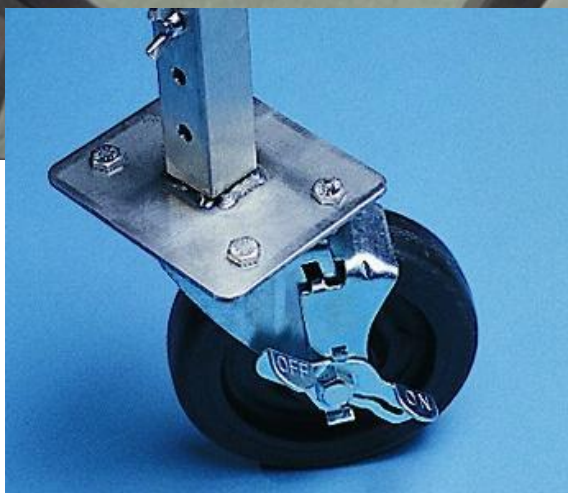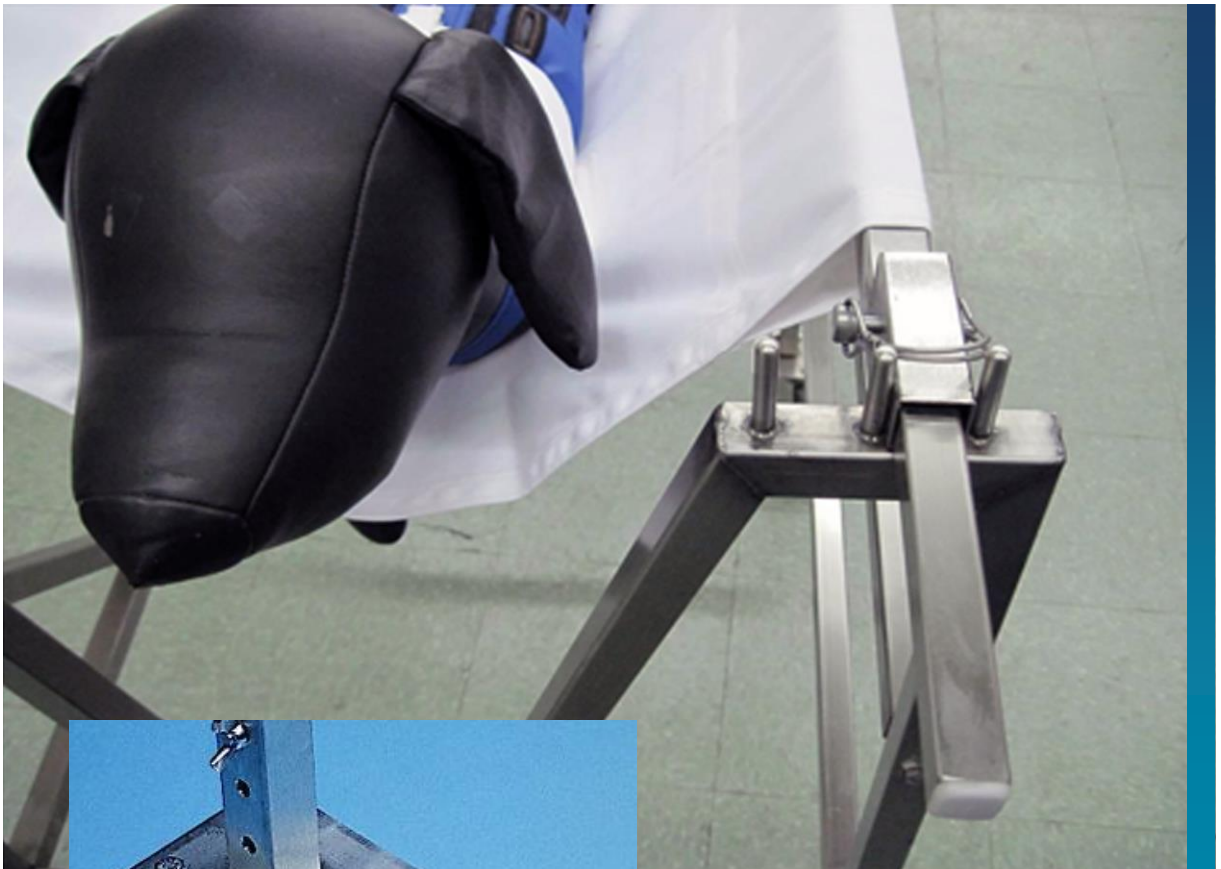

FLAT-TOP FRAMES

FRAME



- For operations using multiple animal at any one time LOMIR manufactures multi-place slings. A flat-top design has the subjects placed into a taut sling cover in either two or four place configurations.

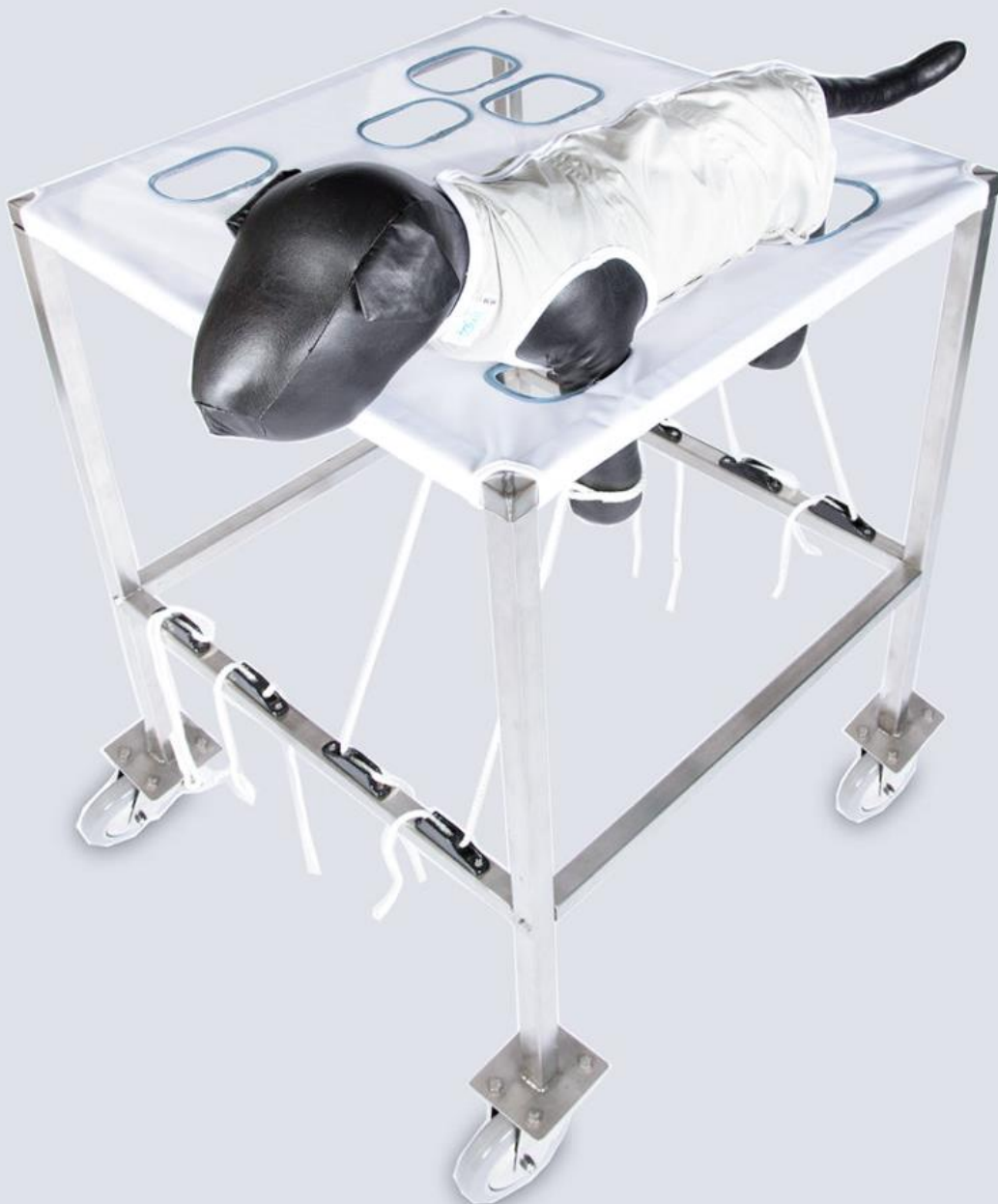

FLAT-TOP FRAMES

SLING COVER


- Interchangeable sling covers are made for individual species.

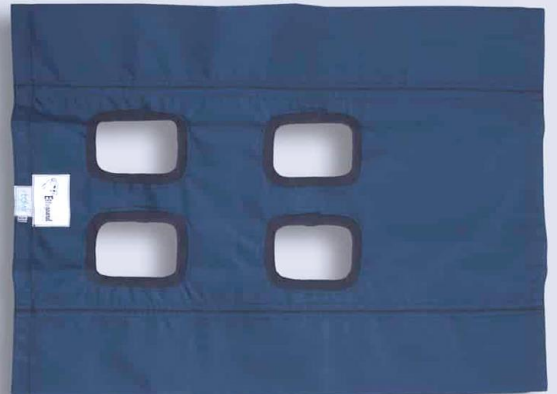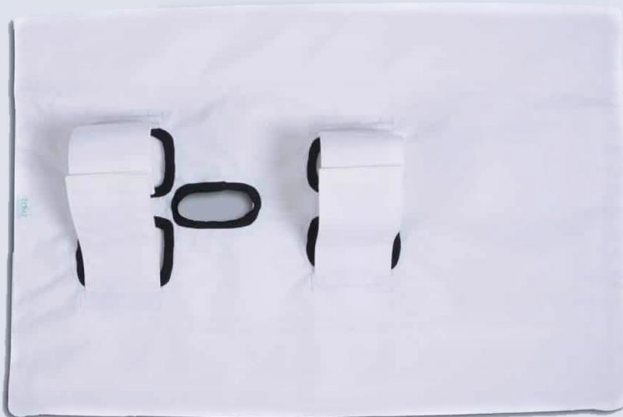

FLAT-TOP FRAMES


SNAP ATTACHMENT

- Cover and frame are matched together using snap fasteners.

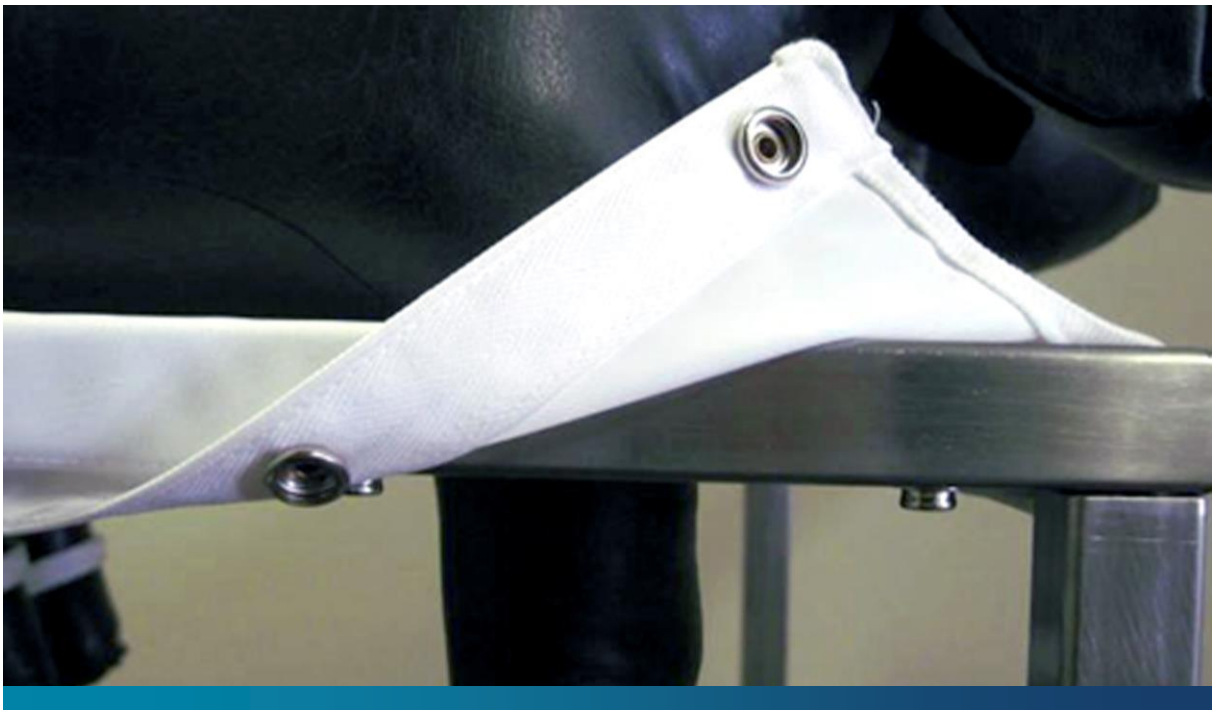

FLAT-TOP FRAMES


FEATURES

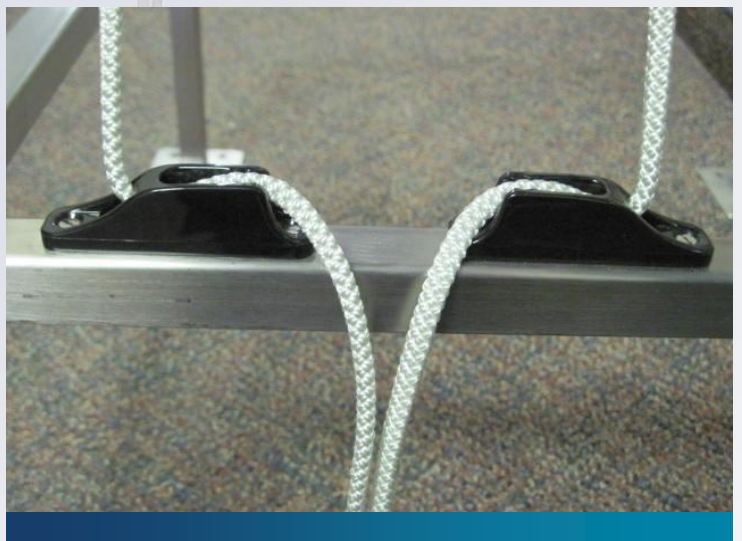

FLAT-TOP FRAMES

FEATURES

- All openings are edged in cotton jersey for comfort.

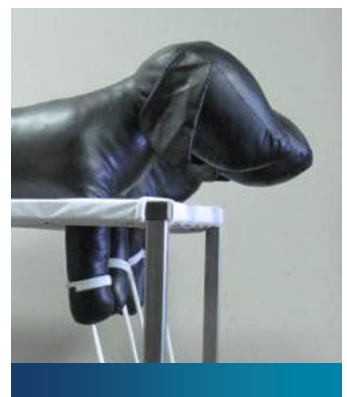

PORTABLE PANEPINTO SLING

OVERVIEW

- Humane, user-friendly, portable restraint
- Developed in partnership with Panepinto Associates
- For use with small to medium sized animals such as:
 - micro/mini pigs
 - Dogs
 - Goats
 - Lambs
- Comes with instructions for:
 - Assembly
 - Storage
 - Handling
- Easy set up and dismantling:
 - No tools required
 - Clevis pins lock the frame in place
- For storage:
 - Fold close and hang on hooks or
 - Use included storage bag



PORTABLE PANEPINTO SLING

FRAME


- To maintain a light weight with maximum strength the frame is constructed from square tubular stainless steel.
- Deep “V” shaped opening at the front and back permit unimpeded access to the animal for easy sampling and instrumentation.



PORTABLE PANEPINTO SLING

SLING COVER

- The sling is machine washable. All openings on the sling are trimmed with a cotton jersey fabric for comfort. Velcro straps closing over the back provide for quick and secure restraint while maintaining adjustable fit for individual animals.

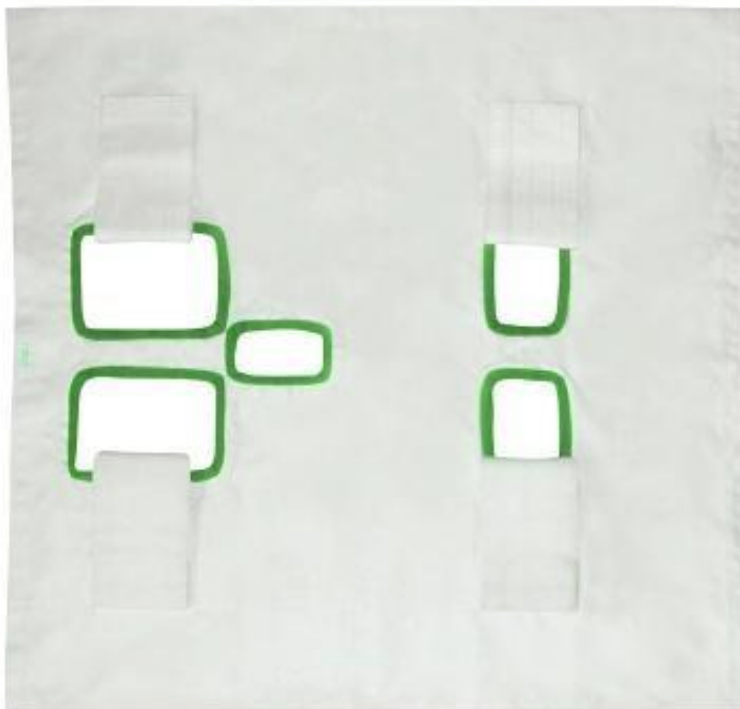



DESIGN. INTEGRITY. RELIABILITY.

**For more information, or to order please visit
us online at www.lomir.com or call 1-877-425-3604**

Lomir Biomedical Inc.

95 Huot, Notre-Dame-de-l'Île-Perrot, QC
Canada J7W 1M3
Tel: 514-425-3604
Toll Free: 1-877-425-3604
Fax: 514-425-3605
Email: info@lomir.com

Lomir Biomedical Inc.

C/O Agenda Life Sciences
UK PO Box 24 Hull
HU128YK, UK
Tel: 03456 44 55 45
Email: info@lomir.com

Lomir Inc.

PO Box 778
Malone, NY
USA 12953
Tel: 518-483-7697
Toll Free: 1-877-425-3604
Email: info@lomir.com