



P.E.T 2028

CARE & USAGE

PRIMATE ENRICHMENT

Thank you for your purchase of the 2028 P.E.T. This enrichment device is designed to be used with New World Primates. The device is manufactured from food grade polypropylene, all fasteners are stainless steel.

CLEANING:

The P.E.T. environmental enrichment device may be cleaned by a number of techniques. For example, emersion in a bleach or disinfectant solution followed by rinsing, placing in a cage or tunnel washer. They may be sterilized using an autoclave.

FILLING:

For filling and cleaning the 2028 P.E.T., unscrew the front and remove the disk. Use a tongue depressor to fill cavities. For sauce, jelly and gels use a large spatula to quickly fill all cavities.

FOOD/TREATS:

The 2028 device can be filled with a range of food items. The following are suggestions: dried fruits, fruit sauce or jellies, nut butter, propriety treats or diet (e.g. Lab Diet's Callitrichid diets 5LK6 or 5LK7). Clear H₂O gels such as LabGel® could also be used.



Should you have any questions or need more information please contact our Customer Service Department toll free in North America at 1-877-425-3604 or send an e-mail to info@lomir.com.

Lomir Biomedical Inc.

95 Huot, Notre-Dame-de-l'Île-Perrot, QC
Canada J7W 1M3
Tel: 514-425-3604
Toll Free: 1-877-425-3604
Fax: 514-425-3605
Email: info@lomir.com

Lomir Biomedical Inc.

C/O Agenda Life Sciences
UK PO Box 24 Hull
HU128YK, UK
Tel: 03456 44 55 45
Email: info@lomir.com

Lomir Inc.

PO Box 778
Malone, NY
USA 12953
Tel: 518-483-7697
Toll Free: 1-877-425-3604
Email: info@lomir.com



P.E.T 2028

CARE & USAGE

PLACEMENT OF DEVICE:

The 2028 can be hung inside an enclosure or cage or can be mounted on the front of a cage. All devices are supplied with two stainless steel eye bolts for mounting, an additional set of two threaded holes at the base of the device allow for additional attachments.

To dismantle for cleaning, changing the plates and filling with food/treats, hold stainless hex bolt with a wrench and unscrew plastic disk on the outer portion of the device.

Animals may use various tools to forage in the smaller holes located in the top and bottom trays and when the plate with the smaller holes is used.



Should you have any questions or need more information please contact our Customer Service Department toll free in North America at 1-877-425-3604 or send an e-mail to info@lomir.com.

Lomir Biomedical Inc.

95 Huot, Notre-Dame-de-l'Île-Perrot, QC
Canada J7W 1M3
Tel: 514-425-3604
Toll Free: 1-877-425-3604
Fax: 514-425-3605
Email: info@lomir.com

Lomir Biomedical Inc.

C/O Agenda Life Sciences
UK PO Box 24 Hull
HU128YK, UK
Tel: 03456 44 55 45
Email: info@lomir.com

Lomir Inc.

PO Box 778
Malone, NY
USA 12953
Tel: 518-483-7697
Toll Free: 1-877-425-3604
Email: info@lomir.com