



P.E.T 2040

CARE & USAGE

PRIMATE ENRICHMENT

Thank you for your purchase of the 2040 P.E.T. This enrichment device is designed to be used with New World Primates. The device is manufactured from food grade polypropylene, all fasteners are stainless steel.

CLEANING:

The P.E.T. environmental enrichment device may be cleaned by a number of techniques. For example, emersion in a bleach or disinfectant solution followed by rinsing, placing in a cage or tunnel washer. They may be sterilized using an autoclave.

FILLING:

Use a bag to fill individual cavities.

FOOD/TREATS:

The 2040 device can be filled with a range of food items. The following are suggestions: dried fruits, fruit sauce or jellies, nut butter, propriety treats or diet (e.g. Lab Diet's Callitrichid diets 5LK6 or 5LK7). Clear H2O gels such as LabGel® could also be used.

PLACEMENT OF DEVICE:

The 2040 can be hung inside an enclosure or cage or can be mounted on the front of a cage. All devices are supplied with two stainless steel eye bolts for mounting, an additional set of two threaded holes at the base of the device allow for additional attachments.



Should you have any questions or need more information please contact our Customer Service Department toll free in North America at 1-877-425-3604 or send an e-mail to info@lomir.com.

Lomir Biomedical Inc.

95 Huot, Notre-Dame-de-l'Île-Perrot, QC
Canada J7W 1M3
Tel: 514-425-3604
Toll Free: 1-877-425-3604
Fax: 514-425-3605
Email: info@lomir.com

Lomir Biomedical Inc.

C/O Agenda Life Sciences
UK PO Box 24 Hull
HU128YK, UK
Tel: 03456 44 55 45
Email: info@lomir.com

Lomir Inc.

PO Box 778
Malone, NY
USA 12953
Tel: 518-483-7697
Toll Free: 1-877-425-3604
Email: info@lomir.com