



CARE INSTRUCTIONS

FOR YOUR SNUGGLES

To clean SNUGGLES, start by soaking for up to an hour in a quaternary ammonium disinfectant, this removes heavy soiling, any odor as well as disinfecting. Do not use chlorine bleach. Then proceed to wash using industrial detergent (such as cage wash detergent) or laundry soap in a washing machine.

Similar results can be achieved by placing SNUGGLES in a cage washer. All disinfectants and detergents are used at the manufacturers' recommended concentrations.

Temperature – Best results will be achieved in warm water.

For best results, close any Velcro fasteners, this protects the Velcro from pulling or attaching to other fabrics and SNUGGLES. SNUGGLES may be machine dried.

All SNUGGLES are protected by US patent # 5,975,028. All rights are reserved.



Should you have any questions or need more information please contact our Customer Service Department toll free in North America at 1-877-425-3604 or send an e-mail to info@lomir.com.

Lomir Biomedical Inc.

95 Huot, Notre-Dame-de-l'Île-Perrot, QC
Canada J7W 1M3
Tel: 514-425-3604
Toll Free: 1-877-425-3604
Fax: 514-425-3605
Email: info@lomir.com

Lomir Biomedical Inc.

C/O Agenda Life Sciences
UK PO Box 24 Hull
HU128YK, UK
Tel: 03456 44 55 45
Email: info@lomir.com

Lomir Inc.

PO Box 778
Malone, NY
USA 12953
Tel: 518-483-7697
Toll Free: 1-877-425-3604
Email: info@lomir.com